

# Important notes

## The return transport - one call is enough!

The return transport is carried out by our logistics partner.

Please observe the export regulations of your country.

GO! Express & Logistics GmbH

Customer-Nr: 103977

Phone Germany: 0800 1112828

Phone abroad: +49 7131 155650

## Please note on the day of sampling:

- Registration of the collection until 12 o'clock at the latest
- Collection takes place until 2 pm
- Last possible return day is always Thursday
- No collection before holidays

## Content:

- 20 Rodac plates for the examination of surfaces
- 8 Bioindicators in blister packaging
- 8 empty sterile tubes
- 1 Transport control bioindicator
- 1 sterile tweezers
- 1 waterproof pen
- 1 Bottle of hand sanitizer
- 1 Cooling element

## Do you have any questions:

Write to us at:  
[info@hymo-box.de](mailto:info@hymo-box.de)




For more information,  
please visit:  
[www.hymo-box.de](http://www.hymo-box.de)

Kitchen Check

# Hygiene monitoring made easy

The  
HyMo-  
Box



Take your hygiene management into your own hands with the HyMo-Box!

The HyMo box contains bioindicators that you simply add during the dishwashing process. These provide them with reliable findings on the cleaning and disinfection performance of your dishwasher.

In addition, surface testing can be used to determine the Hygiene can be determined directly on the wash ware, but also on surfaces in the immediate vicinity.

Please note the instructions for return transport on the reverse side.

# Step by step guide

## for the HyMo box kitchen check

Your benefits: You can easily and reliably monitor your cleaning and disinfection measures. It's all inclusive - no hidden costs! Our DAkkS accreditation guarantees you the highest quality of our services.



1

- Open HyMo-Box.
- Place cooling element in the refrigerator.
- In case of later processing, store the box with the RODAC plates and the bioindicators also in the refrigerator.



3

- Disinfect hands with the enclosed Disinfect hands with disinfectant.

2

- Define day of performance of tests and organize return transport for the same day.
- Attention: registration of the return transport must be made before 12 noon, the collection of the shipment will be made until 2 pm, no return transport on Fridays.




4

- Remove the 8 bioindicators from the packaging. To do this, tear open the blister packaging at the top with your hands, hold the bottom end of the packaging with one hand and let the bioindicator slide into the compartments of the cutlery basket without touching it.
- It is also possible to place the bioindicator between plates or in other relevant places in the rinsing chamber of the machine. Ensure that it is adequately fastened.
- Transport control remains untouched in the box.



5

- Start the machine with the program to be checked.
- Disinfect hands again.
- At the end of the program, transfer the bio-indicators individually into an empty test tube using tweezers and close tightly.
- When taking samples from the wash ware, touch it only at the edge. Allow to cool.

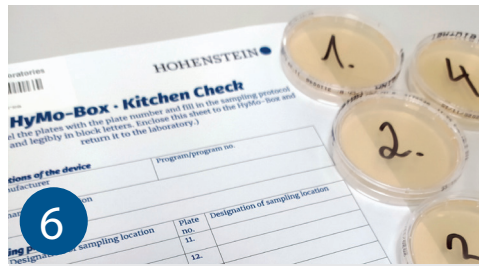

6

#### PROBATE:

- Define sampling sites for surface examination.
- Select items to be washed (ideally 10 different pieces of crockery and cutlery) from different places in the dishwasher.
- Furthermore, the interior of the machine as well as sensitive areas of the environment can be selected.
- Enter these in the sampling plan. Label the RODAC plates with numbers according to the sampling plan.



7

- Disinfect hands.
- Open the lid of a RODAC plate.
- Press the culture medium surface for approx. Press lightly on the area to be examined for approx. 5 seconds - do not move the RODAC plate if possible.
- Close the lid again.
- With all further RODAC plates proceed according to your sampling plan.



8

- Stack the RODAC plates back into the box.
- Put the box, the 8 tubes with the bioindicators and the transport control as well as the cooling element back into the box and close it.
- The sampling plan is placed on the box and the outer box is closed.