

Important notes

The return transport - one call is enough!

The return transport is carried out by our logistics partner.

Please observe the export regulations of your country.

GO! Express & Logistics GmbH

Customer-Nr: 103977

Phone Germany: 0800 1112828

Phone abroad: +49 7131 155650

Please note on the day of sampling:

- Registration of the collection until 12 o'clock at the latest
- Collection takes place until 2 pm
- Last possible return day is always Thursday
- No collection before holidays

Content:

- RODAC PLATES TO SURFACE
- INVESTIGATIONS:
 - Surface check: 20 pieces or
 - Surface check XL: 40 pieces
- 1 bottle of hand sanitizer
- 1 Cooling element
- 1 waterproof pen

Do you have any questions:

Write to us at:
info@hymo-box.de



For more information,
please visit:
www.hymo-box.de

Surface Check & Surface Check XL

Hygiene monitoring made easy

The
HyMo-
Box



Take your hygiene management into your own hands with the HyMo-Box!

The HyMo-Box provides you with reliable findings, whether your focus is on the hygienic safety of work surfaces, floors or fittings.

Using the globally established smear test method, our accredited microbiological laboratory determines important key figures and recommendations for all surfaces examined.

Please note the instructions for return transport on the reverse side.

Step by step guide

for the HyMo surface box

Your advantages: You can easily and reliably monitor your cleaning and disinfection measures. It's all inclusive - there are no hidden costs for you. Our DAkkS accreditation guarantees you the highest quality of our services.



1

- Open HyMo-Box.
- Place cooling element in the refrigerator.
- You can also store the box of RODAC plates in the refrigerator for later sampling.



2

- Define the day of your sampling and organize return transport for the same day accordingly. Attention: Registration of the return transport must be done before 12 noon, the collection of the shipment will be done until 2 pm, no return transport on Fridays.
- Determine sampling locations and enter them in the attached sampling protocol record.



3

- Label the bottom of each RODAC plate according to the sampling plan with the corresponding numbers.



4

- Disinfect hands with the enclosed Disinfect hands with disinfectant.



5

SAMPLING:

- Open the lid of a RODAC plate.
- Press the culture medium surface for approx. 5 seconds under slight pressure on the area to be examined - do not move the RODAC plate if possible.
- Close the lid again.



6

- Proceed with all further RODAC plates according to your sampling plan.
- Stack the RODAC plates back into the box.



7

- Return the sealed box to the box.
- Add cooling element and close box.



8

- Sampling protocol on the box and tape the glue shut.